

# nXDS Pump Accessories

INSTRUCTION MANUAL

A73501840\_D Original instructions

### **Copyright notice**

©Edwards Limited 2024. All rights reserved.

Published: 2/6/2024

### **Associated publications**

#### **Publication title**

**Publication number** 

Vacuum pump and vacuum system safety

P400-40-100

### **Trademark credit**

Edwards and the Edwards logo are trademarks of Edwards Limited, Innovation Drive, Burgess Hill, West Sussex, UK, RH15 9TW

### **Disclaimer**

The content of this manual may change from time to time without notice. We accept no liability for any errors that may appear in this manual nor do we make any expressed or implied warranties regarding the content. As far as practical we have ensured that the products have been designed and constructed to be safe and without risks when properly installed and used in accordance with their operating instructions.

We accept no liability for loss of profit, loss of market or any other indirect or consequential loss whatsoever.

Product warranty and limit of liability are dealt with in our standard terms and conditions of sale or negotiated contract under which this document is supplied.

You must use this product as described in this manual. Read the manual before you install, operate, or maintain the product.

### **Contents**

1.	Safety and compliance	7
	1.1 Definition of Warnings and Cautions	7
	1.2 Trained personnel	7
	1.3 Safety symbols	8
2.	Introduction	9
	2.1 Scope and definitions	. 9
	2.2 Installation and operation safety	9
	2.3 Unpack and inspect	9
	2.4 Disposal of accessories	
	2.5 Spares and accessories	10
3.	Gas ballast adaptor blank	12
	3.1 Description	12
	3.2 Installation	12
	3.2.1 Install the gas ballast adaptor blank	12
4.	Gas ballast adaptor	13
	4.1 Description	13
	4.2 Technical data	13
	4.3 Installation	13
	4.3.1 Install the gas ballast adaptor	14
5.	Inlet/exhaust filter	16
	5.1 Description	16
	5.2 Technical data	16
	5.3 Installation	17
	5.3.1 Install the inlet/exhaust filter	17
	5.4 Maintenance	17
	5.4.1 Introduction	
	5.4.2 Removal of elements	17
	5.4.3 Inspection and replacement of the element	18
6.	Silencer	19
	6.1 Description	19
	6.1.1 Application	19
	6.1.2 Connections	19
	6.2 Technical data	19
	6.3 Installation	20
	6.3.1 Install the nXDS silencer	20

	6.4 Maintenance	21
	6.4.1 Introduction	21
	6.4.2 Removal of elements	21
	6.4.3 Inspection of elements	21
	6.4.4 Replacement of elements	22
<b>7.</b>	Exhaust nozzle	23
	7.1 Description	23
	7.2 Technical data	23
	7.3 Installation	23
	7.3.1 Install the exhaust nozzle	23
8.	Mountable vibration isolators	
	8.1 Description	
	8.2 Technical data	25
	8.3 Installation	
	8.3.1 Install the Edwards mountable vibration isolators	26
9	Chemical resistance conversion kit	27
٠.	9.1 Description	
	9.2 Installation.	
	9.2.1 Install the chemical resistance conversion kit	
	9.2.2 Inlet/exhaust port conversion	
	3.2.2 Interpexhaust port conversion	23
10	. Pump-to-controller cable	30
	10.1 Description	30
	10.2 Technical data	
	10.3 Installation	
		_
11	. Legal declarations	31

## **List of Figures**

Figure 1: Accessories for nXDS pumps - general view (o-ring seals not shown)	11
Figure 2: Gas ballast adaptor blank fitting	12
Figure 3: Gas ballast adaptor conversion	14
Figure 4: Inlet/exhaust dust filter components and dimensions	16
Figure 5: Silencer dimensions	20
Figure 6: Nozzle and dimensions	23
Figure 7: Vibration isolator dimensions	25
Figure 8: Inlet/exhaust port conversion	29

### **List of Tables**

Table 1: Gas ballast adaptor blank - checklist of components (A735-01-806)	12
Table 2: Gas ballast adaptor technical data	13
Table 3: Gas ballast adaptor - checklist of components	13
Table 4: Inlet/exhaust filter technical data	16
Table 5: Table: Inlet/exhaust filter - checklist of components NW25 (A505-97-805)/ NW40 (A505-97-806)	١
	17
Table 6: NW25/NW40 filter element replacement kit	18
Table 7: Silencer technical data	19
Table 8: Silencer - checklist of components (A505-97-000)	20
Table 9: Silencer element replacement kit	22
Table 10: Nozzle technical data	23
Table 11: Exhaust nozzle - checklist of components (A505-09-000)	23
Table 12: Vibration isolator technical data	25
Table 13: Mountable vibration isolators - checklist of components (A248-01-441)	26
Table 14: Chemical resistance conversion kit - checklist of components nXDS6i/10i/15i (A735-01-807).	27
Table 15: Chemical resistance conversion kit - checklist of components nXDS20i (A735-01-808)	27
Table 16: Ordering information	30
Table 17: Checklist of components	30

### 1. Safety and compliance

For safe operation from the start, read these instructions carefully before you install or commission the equipment and keep them safe for future use. Read all the safety instructions in this section and the rest of this manual carefully and make sure that you obey these instructions.

The instruction manual is an important safety document that we often deliver digitally. It is your responsibility to keep the instruction manual available and visible while working with the equipment. Please download the digital version of the instruction manual for use on your device or print it if a device will not be available.

### 1.1 Definition of Warnings and Cautions

Important safety information is highlighted as warning and caution instructions which are defined as follows. Different symbols are used according to the type of hazard.

#### **WARNING:**

If you do not obey a warning, there is a risk of injury or death.

#### **CAUTION:**

If you do not obey a caution, there is a risk of minor injury, damage to equipment, related equipment or process.

#### **NOTICE:**

Information about properties or instructions for an action which, if ignored, will cause damage to the equipment.

We reserve the right to change the design and the stated data. The illustrations are not binding.

### 1.2 Trained personnel

For the operation of this equipment "trained personnel" are:

- skilled workers with knowledge in the fields of mechanics, electrical engineering, pollution abatement and vacuum technology and
- personnel specially trained for the operation of vacuum pumps

### 1.3 Safety symbols

The safety symbols on the products show the areas where care and attention is necessary.

The safety symbols that we use on the product or in the product documentation have the following meanings:



#### Warning/Caution

Risk of injury and/or damage to equipment. An appropriate safety instruction must be followed or a potential hazard exists.



#### Warning - Use protective equipment

Risk of injury. Use appropriate Personal Protective Equipment (PPE) when performing the task.

### 2. Introduction

#### 2.1 Scope and definitions

This manual provides installation, operation and maintenance instructions for the Edwards range of accessories for the nXDS pump range.

The accessories are shown in *Figure: Accessories for nXDS pumps - general view (o-ring seals not shown)* on page 11. The Item Numbers for the accessories are listed in the appropriate sections. The accessories must be used as specified in this manual. Read this manual before installing accessories onto the nXDS pump.

#### 2.2 Installation and operation safety



#### **WARNING: INSTALLATION SAFETY**

Risk of injury or damage to the equipment. Use the procedures described in this manual to install the accessory. Obey all safety instructions and take note of all appropriate precautions. If not, damage to the accessory or other equipment and injury to people can result.



#### **WARNING: OPERATION SAFETY**

Risk of injury or damage to the equipment. The user of the nXDS pump system is responsible for the safe operation and monitoring of the system.



#### **WARNING: INSTALLATION SAFETY**

Risk of injury or damage to the equipment. Before installing the accessory, ensure the pump is off and the controller is isolated as described below.

Before installing the accessory:

- Switch off the nXDS pump and wait until the pump has stopped rotating.
- Isolate the pump controller from the electrical supply.

### 2.3 Unpack and inspect

Remove all packing materials and check the accessory. If it is damaged, notify the supplier and the carrier in writing within three working days; state the item number of the accessory together with the order number and the suppliers invoice number. Retain all packing materials for inspection. Do not use the accessory if it is damaged.

Check that the packaging contains the items shown in the Checklist of items table. If any of the items are missing, notify the supplier in writing within three working days.

### 2.4 Disposal of accessories

Dispose of the used accessories and accessory spares safely in accordance with all local and national safety environmental requirements.


### 2.5 Spares and accessories

Edwards products, spares and accessories are available from Edwards companies in Belgium, Brazil, Canada, France, Germany, Hong Kong, Italy, Japan, Korea, Switzerland, United Kingdom, U.S.A, and a worldwide network of distributors. The majority of these centers employ service Engineers who have undergone comprehensive Edwards training courses.

Order spare parts and accessories from the nearest Edwards company or distributor. When ordering, please state for each part required:

- Model and Item Number of the equipment
- Serial number (if any)
- Item Number and description of part.

Figure 1 Accessories for nXDS pumps - general view (o-ring seals not shown)

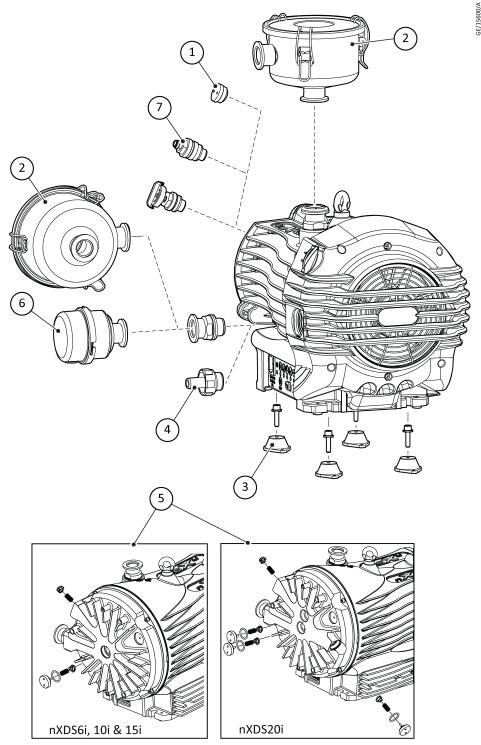

- 1. Gas ballast adaptor blank
- 3. Vibration isolators
- 5. Chemical resistance conversion kit
- 7. Gas ballast adaptor

- 2. Inlet/exhaust filter
- 4. Exhaust nozzle
- 6. Silencer
- 8. Pump to controller cable (not shown)

### 3. Gas ballast adaptor blank

#### 3.1 Description

The gas ballast adaptor blank (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown)* on page 11, item 1) allows conversion from a standard nXDS pump into an nXDS-R pump. The gas ballast control is replaced with a blank adaptor to prevent accidental air admittance into the pump. This feature is also useful for applications such as rare gas recirculation or gas recovery.

#### 3.2 Installation

Table 1 Gas ballast adaptor blank - checklist of components (A735-01-806)


Qty	Description	Check (✓)
1	Gas ballast adaptor blank	
1	O-ring 19.6ID X 2.4	

#### 3.2.1 Install the gas ballast adaptor blank

The method for installing the gas ballast adaptor blank follows some of the stages used to install the gas ballast adaptor. To install the gas ballast adaptor blank, refer to stages 1, 2, 3, 4 and 6 of the gas ballast adaptor fitting instructions in *Install the gas ballast adaptor* on page 14. For the instructions on how to fit the adaptor blank refer to the instruction and *Figure: Gas ballast adaptor blank fitting* on page 12 below.

Using a pin spanner (not provided), fit the adaptor blank and tighten to 6 Nm.

Figure 2 Gas ballast adaptor blank fitting

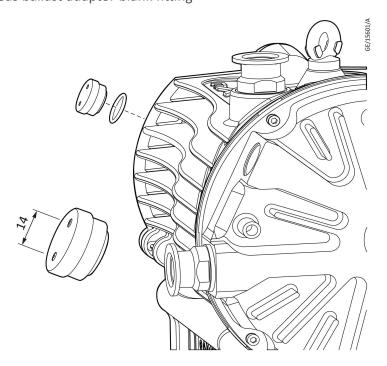

### 4. Gas ballast adaptor

### 4.1 Description

The gas ballast adaptor (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown*) on page 11, item 7) fits in place of the gas ballast control. The gas ballast adaptor allows connection of a controlled supply of inert gas such as nitrogen into the low vacuum stage of the pump.

The gas purge can be used to dilute flammable gases down to a safe level (Edwards recommends 1/4 of the Lower Explosive Limit). An inert purge can also be used to reduce the likelihood of gas vapours such as iodine from condensing at the outlet stage of the pump.

#### 4.2 Technical data

Table 2 Gas ballast adaptor technical data

Part number	Description	Maximum supply pressure	GB flow
A73501809	A73501809 Gas ballast adaptor with fine restrictor		0.5 l/m
A73501811 Gas ballast adaptor without fine restrictor		1 bara	Standard*

<sup>\*</sup> Refer to the n instruction manual for standard flow rate.


#### 4.3 Installation

Table 3 Gas ballast adaptor - checklist of components


Qty	Description	Check (√)
1	Gas ballast adaptor	
1	O-ring 19.6 ID x 2.4	


#### 4.3.1 Install the gas ballast adaptor

Figure 3 Gas ballast adaptor conversion

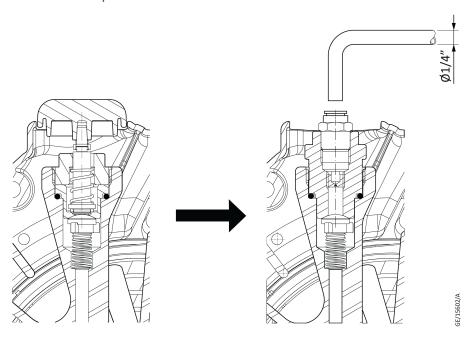

Ensure the pump is switched off before fitting.

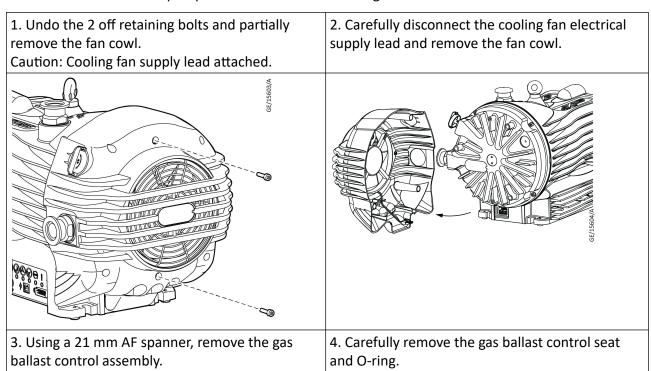

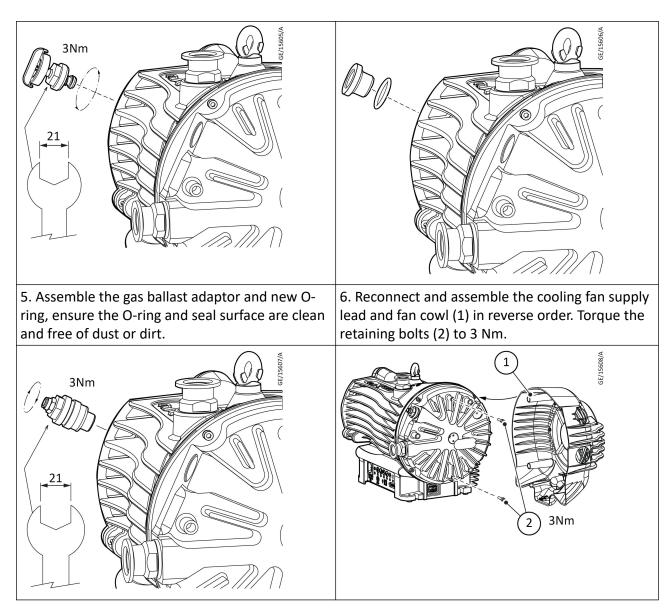

As a precaution, Edwards recommends that a complete a vacuum leak check be performed after this installation.

#### Tip:

- Ensure the cooling fan operates correctly following the installation of the gas ballast adaptor.
- Retain the gas ballast control assembly and seat for future use if required.

### 5. Inlet/exhaust filter

#### 5.1 Description

#### **CAUTION: OPERATION SAFETY**



Risk of damage to the equipment. This filter is not designed to withstand positive pressures. Ensure that any system valves are correctly scheduled to avoid pumping to a closed exhaust line. Refer to the Installation section of the nXDS pump instruction manual.


The inlet/exhaust filter performs a number of functions. It can be connected in-line between the nXDS pump exhaust and the exhaust extraction system (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown)* on page 11, item 2), this is especially useful for process applications such as gas recirculation. Alternatively, it can be connected in-line between the nXDS pump inlet and the chamber outlet connection (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown)* on page 11, item 2) to prevent particles from migrating (back streaming) into the system, or to stop particulate from entering the pump. The filter is supplied ready to use with a 5 micron element. A 1 micron element can be purchased separately and used to replace the 5 micron element if required.

#### 5.2 Technical data

Table 4 Inlet/exhaust filter technical data

Inlet/outlet port size	NW25 (A505-97-805)
	NW40 (A505-97-806)
Leak tightness	1 x 10-5 mbar.ls-1

Figure 4 Inlet/exhaust dust filter components and dimensions

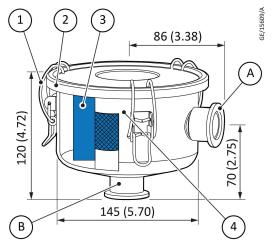

- A. Inlet
- 1. Lid fastening clamp
- 3. Polyester element

- B. Exhaust
- 2. Filter lid
- 4. Filter body

#### 5.3 Installation

Table 5 Table: Inlet/exhaust filter - checklist of components NW25 (A505-97-805)/ NW40 (A505-97-806)

Qty	Description	Check (✓)
1	Filter assembly	
2	NW clamp	
2	Centring ring	

#### 5.3.1 Install the inlet/exhaust filter

Ensure the pump is switched off before fitting.

- 1. Inspect the mating port and seals faces, ensure they are free of dirt or any particulate which could cause the vacuum seal to leak.
- 2. Place the centering ring in the end of the NW25 filter port and seat the filter onto the pump mating port. Ensure the filter is fitted in the correct orientation (refer to Figure: Accessories for nXDS pumps general view (o-ring seals not shown) on page 11 and Figure: Inlet/exhaust dust filter components and dimensions on page 16).
- 3. Fit the NW25 retaining clamp and tighten to secure the filter.

#### 5.4 Maintenance

#### 5.4.1 Introduction

Edwards recommends that a filter element replacement (*Figure: Inlet/exhaust dust filter components and dimensions* on page 16, item 3) should coincide with a tip seal change. Refer to the maintenance plan in the nXDS pump manual for recommendations on changing the tip seal.

#### **5.4.2** Removal of elements



#### **WARNING: USE PROTECTIVE EQUIPMENT**

Risk of injury or damage to the equipment. Take all necessary precautions if toxic or dangerous substances have been pumped. Wear PPE when handling contaminated filters and filter elements.

#### WARNING: OPERATION SAFETY



Risk of injury or damage to the equipment. Do not attempt to dismantle the filter whilst the pump is running. Ensure the pump is isolated from its electrical supply before removing the filter.

Do not inhale the tip seal dust. Do not blow the tip seal dust from the filter with compressed air.

Remove the filter from the system. Release the clamps (Figure: Inlet/exhaust dust filter components and dimensions on page 16, item 1) and remove the lid (Figure: Inlet/

exhaust dust filter components and dimensions on page 16, item 3) from the filter body (Figure: Inlet/exhaust dust filter components and dimensions on page 16, item 4). Remove the element for inspection.

#### 5.4.3 Inspection and replacement of the element



#### **WARNING: USE PROTECTIVE EQUIPMENT**

Risk of injury or damage to the equipment. Take all necessary precautions if toxic or dangerous substances have been pumped. Wear PPE when handling contaminated filters and filter elements.



#### **WARNING: HAZARDS GASES**

Risk of injury or damage to the equipment. At high temperature (in excess of 250 °C) tip seal wear product will begin to decompose, giving rise to gaseous fumes which can produce unpleasant symptoms.

The filter may be contaminated with the process chemicals that have been pumped during operation. Ensure that the pump is decontaminated before maintenance and that adequate precautions are taken to protect people from the effects of dangerous substances if contamination has occurred.

Renewal of the element is recommended regardless of the level of contamination. If the element is heavily contaminated with tip-seal dust it must be replaced or cleaned in order to avoid a build up of dust that may adversely affect the pump performance. Rinse the element if necessary with lukewarm water and mild detergent. Observe local guidelines to advise the safe disposal of PTFE dust.

Fitting the new element is the reverse of removal.

Table 6 NW25/NW40 filter element replacement kit

Product description	Ordering information
5 micron element	A505-97-802
1 micron element	A505-97-803

#### 6. Silencer

#### 6.1 Description

The nXDS silencer has two main functions.

- To reduce exhaust noise during pump down and gas ballast operation.
- To collect and retain tip seal dust as the pump is operated.

Exhaust gases exit the exhaust port when the pump is being used to evacuate a process chamber, when the gas ballast control is being used, or both. The noise level when pumping down rises as the throughput increases and the silencer is used to reduce this noise level.

The outer can is made of pressed steel and is split to allow access to the filter elements.

The filter element housing is made from Santoprene® thermoplastic elastomer. The element itself is polyester. The foam block is manufactured from filter foam.

#### ■ Note:

These elements can only be used in dry applications and must not be allowed to moisten.

A pressure relief valve is not fitted to the silencer.

#### 6.1.1 Application



#### **WARNING: INSTALLATION SAFETY**

Risk of injury or damage to the equipment. Our nXDS silencer is not recommended for use with pumping hazardous substances.

#### 6.1.2 Connections


The nXDS silencer has been manufactured with an NW25 fitting to enable fitting to the nXDS pump exhaust (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown*) on page 11, item 6). An NW25 clamp and centering ring are included. A sectional view of the silencer is shown in *Figure: Silencer dimensions* on page 20.

#### 6.2 Technical data

Table 7 Silencer technical data

Maximum back pressure	0.2 bar gauge
Tip seal dust	Pump performance and silencer performance is unaffected providing elements are serviced in accordance with recommendations.
Dimensions	Refer to Figure: Silencer dimensions on page 20
Mass	265 g
Port size	NW25

Figure 5 Silencer dimensions

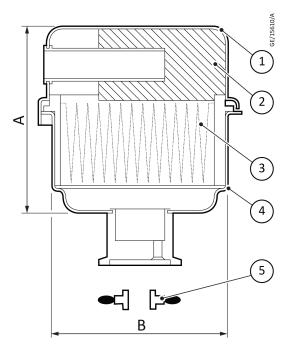

- A. 94 mm
- 1. Top casting
- 3. Polyester element
- 5. Centering ring

- B. 81 mm
- 2. Foam block
- 4. Bottom casing

#### 6.3 Installation

*Table 8 Silencer - checklist of components (A505-97-000)* 

Qty	Description	Check (√)
1	Silencer body	
1	NW25 clamp	
1	Centering ring	

#### 6.3.1 Install the nXDS silencer

Ensure the pump is switched off before fitting the silencer.

- 1. Clean the face of the pump exhaust.
- 2. Place the centering O-ring (*Figure: Silencer dimensions* on page 20, item 5) in the end of the NW25 backing port and seat the silencer in the pump exhaust.
- 3. Fit the clamp and tighten to secure the silencer.

#### 6.4 Maintenance

#### 6.4.1 Introduction

Figure: Inlet/exhaust dust filter components and dimensions on page 16 shows the foam block (item 2) and polyester element (item 3) which need to be inspected periodically. Edwards recommends both are replaced every 9000 hours of pump use. The silencer has been designed to handle 10 cc of tip-seal dust without affecting the performance of the pump. Replacement kit A505-97-800 contains both items.

#### 6.4.2 Removal of elements



#### **WARNING: USE PROTECTIVE EQUIPMENT**

Risk of injury or damage to the equipment. Take all necessary precautions if toxic or dangerous substances have been pumped. Wear PPE when handling contaminated filters and filter elements.





Risk of injury or damage to the equipment. Do not attempt to dismantle the silencer whilst the pump is running. Ensure the pump is isolated from its electrical supply before removing the silencer.

Do not inhale the tip seal dust. Do not blow the tip seal dust from the silencer with compressed air.

Ensure the pump is switched off before removing the silencer.

Remove the NW25 clamp and silencer from the pump exhaust. Hold the top half of the silencer with one hand, and the bottom half with the other hand. Keeping one hand steady, twist clockwise with the other, and the two halves will separate. The polyester element will be retained in the top half. With the element uppermost, lift the element out by lightly gripping the protruding surface of the element. Put the element to one side ready to inspect it. The foam block is pushed onto the outlet pipe and is trapped into position by the polyester element.

#### 6.4.3 Inspection of elements



#### WARNING: USE PROTECTIVE EQUIPMENT

Risk of injury or damage to the equipment. Take all necessary precautions if toxic or dangerous substances have been pumped. Wear PPE when handling contaminated filters and filter elements.



#### **WARNING: HAZARDS GASES**

Risk of injury or damage to the equipment. The tip seal wear product will begin to decompose if the temperature exceeds 250 °C, giving rise to gaseous fumes that can produce unpleasant symptoms.

The silencer will be contaminated with the process chemicals that have been pumped during operation. Ensure that the pump is decontaminated before maintenance and that adequate precautions are taken to protect people from the effects of dangerous substances if contamination has occurred.

Renewal of the elements is recommended regardless of their level of contamination. Neither element should be heavily contaminated with tip seal dust, but if either is, they MUST be replaced or cleaned in order to avoid a build up of dust that may adversely affect performance. Rinse the elements through with soap and water to clean them. Observe local guidelines to advise the safe disposal of PTFE dust.

#### 6.4.4 Replacement of elements

Fitting the new elements is the reverse of removal, fit the foam block first and then the polyester element. Fit both halves of the silencer together and twist anticlockwise to secure. Ensure fastening is complete by looking at the bayonet fitting and checking it is home.

Table 9 Silencer element replacement kit

Product description	Ordering information
Silencer spares kit	A505-97-800

### 7. Exhaust nozzle

#### 7.1 Description


The exhaust nozzle (*Figure: Accessories for nXDS pumps - general view (o-ring seals not shown)* on page 11, item 4) can be used to connect the pump exhaust using a plastic hose in place of a steel bellows. The Nozzle will accept a 12 to 15 mm plastic hose.

#### 7.2 Technical data

Table 10 Nozzle technical data

Nozzle thread	3/4" BSP
Nozzle tube Ø	15 mm

Figure 6 Nozzle and dimensions

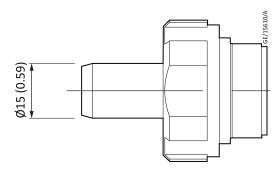

#### 7.3 Installation

*Table 11 Exhaust nozzle - checklist of components (A505-09-000)* 

Qty	Description	Check (√)
1	Nozzle	
1	O-ring	

#### 7.3.1 Install the exhaust nozzle

#### **CAUTION: OPERATION SAFETY**



Risk of damage to the equipment. This exhaust nozzle is not intended to withstand positive pressures. Ensure that any system valves are correctly scheduled to avoid pumping to a closed exhaust line. If using a plastic hose, avoid sharp bends which may kink the hose and restrict the exhaust line. Edwards recommends using reinforced plastic hose. Refer to the installation section of the nXDS pump instruction manual.

Ensure the pump is switched off before fitting.

#### A73501840\_D - Exhaust nozzle

- 1. Using a 30 mm AF spanner (not supplied), remove the NW25 exhaust port from the pump (retain for future use).
- 2. Using a new O-ring provided, manually screw in the plastic exhaust nozzle until resistance can be felt. Tighten the nozzle to 5 Nm using a 36 mm AF spanner (not supplied).
- 3. Push on the plastic hose, use a retaining clip to prevent the hose from becoming unattached from the nozzle.

### 8. Mountable vibration isolators

### 8.1 Description

The Edwards mountable vibration isolators can be used to reduce the transmitted vibration from the pump to a mounting surface such as a system frame.


The mountable vibration isolators have fixing holes to allow the pump to be bolted to the floor or system frame.

#### 8.2 Technical data

Table 12 Vibration isolator technical data

Pump model	Pump mass (kg)	Maximum load per isolator (kg)	Nominal deflection (mm)
nXDS6i	26.2	18.0	4
nXDS10i	25.8	18.0	4
nXDS15i	25.2	18.0	4
nXDS20i	25.6	18.0	4

Figure 7 Vibration isolator dimensions

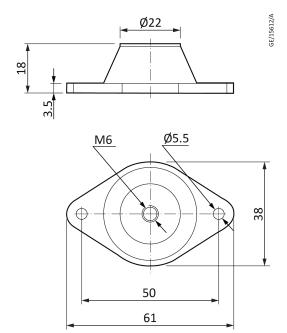

#### 8.3 Installation

Table 13 Mountable vibration isolators - checklist of components (A248-01-441)

Qty	Description	Check (✓)
4	Vibration isolator	
4	Pump fixing screw	
4	Plain washer	
4	Shake-proof washer	

#### 8.3.1 Install the Edwards mountable vibration isolators

Ensure the pump is switched off before fitting.

- 1. Unscrew and remove the standard rubber feet if fitted.
- 2. Fit the vibration isolators to the pump (Figure: Accessories for nXDS pumps general view (o-ring seals not shown) on page 11, item 3) using the screws and washers provided.
- 3. Using 5 mm bolts (not supplied), fix the vibration isolators bolted to the fixing surface or system frame.

Edwards recommends the fitting of flexible tubing or bellows connections between the pump-inlet and the pump-outlet and the rest of the system, to prevent the transmission of stress to the vacuum pipelines and/or equipment.

### 9. Chemical resistance conversion kit

### 9.1 Description

The adaptor kit allows the conversion of a standard nXDS pump into an nXDS-C version. This conversion may be necessary if anticipating the use of the pump on applications involving corrosive substances.

#### 9.2 Installation

Table 14 Chemical resistance conversion kit - checklist of components nXDS6i/10i/15i (A735-01-807)

Qty	Description	Check (✓)
2	Valve pad (Chemraz)	
2	Valve spring	
2	19.6ID O-ring	
2	NW25 stainless steel port	
2	NW25 port O-ring	

Table 15 Chemical resistance conversion kit - checklist of components nXDS20i (A735-01-808)


Qty	Description	Check (✓)
4	Valve pad (Chemraz)	
4	Valve spring	
4	19.6ID O-ring	
2	NW25 stainless steel port	
2	NW25 port O-ring	


#### 9.2.1 Install the chemical resistance conversion kit


Ensure the pump is switched off before fitting. Discard all used 19.6 ID O-ring's. Valve pad conversion

1. Undo the 2 off retaining bolts and partially remove the fan cowl.

Caution: Cooling fan supply lead attached.

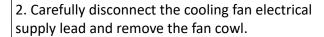


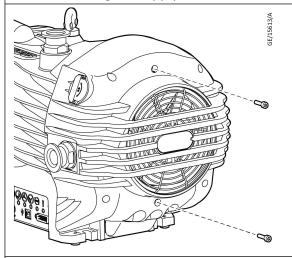


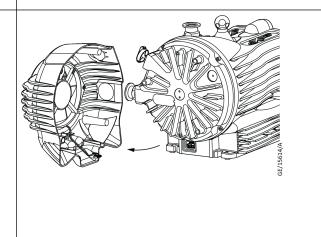


3. Using a 21 mm AF spanner (not supplied), remove the gas ballast control assembly, seat and O-ring.

4a. (nXDS6i,10i &15i) Remove the remaining items (spring and pad) from the gas ballast port. Using a pin spanner, remove the exhaust valve end cap, O-ring, spring and valve pad.

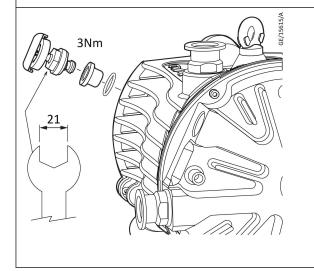

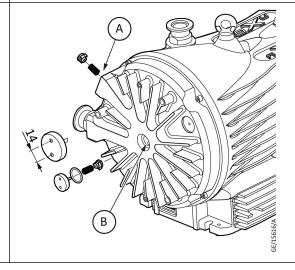

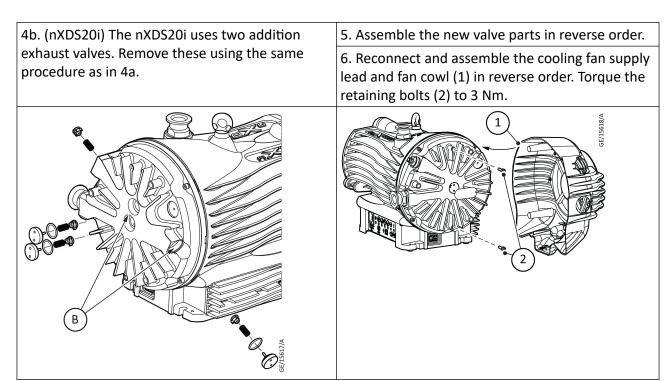


#### A. Gas ballast port

B. Exhaust valves

#### 9.2.2 Inlet/exhaust port conversion

Using a 30 mm AF spanner (not supplied), remove both the inlet and exhaust ports. Locate the stainless steel ports provided in the kit and assemble using new O-rings in reverse order (refer to *Figure: Inlet/exhaust port conversion* on page 29).

Figure 8 Inlet/exhaust port conversion

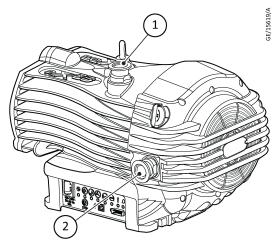

1. Inlet port

2. Exhaust port

### 10. Pump-to-controller cable

### 10.1 Description

The control cable is used to connect the nXDS pump directly to the RS485 or RS232 serial input on the control equipment or a PC, or alternatively, to an Edwards controller such as a TIC (Turbo Instrument Controller) or TAG (Turbo & Active controller).

#### 10.2 Technical data

For full details of the Logic interface technical data, refer to the nXDS pump instruction manual A735-01-880, Section 2.

**Table 16 Ordering information** 

Product description	Ordering information
1 m cable	D397-00-835
2 m cable	D397-00-836
5 m cable	D397-00-837

#### 10.3 Installation

Table 17 Checklist of components

Qty	Description	Check (✓)
1	Control cable	

Ensure the pump is switched off before fitting. Secure the cable using the retaining screws incorporated in the D-type connector.

